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ACCIONES

Aprende a:

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Alertas

TO DO 3

 EQUIPMENT HAS MISSING INFO: ENGINE OF 366215

START

CUSTOMER: ACME Demo

WORKSITE: Houston , TX

UNIT ID: 366215

COMP. DESC.: ENGINE

ASSIGNED BY : Lori_BPNew Beerwart

Add info please

[Comment](#)

 EQUIPMENT HAS MISSING INFO: ENGINE OF 632407

START

CUSTOMER: Acme Demo

WORKSITE: Worksite 2

UNIT ID: 632407

COMP. DESC.: ENGINE

ASSIGNED BY : Lori_BPNew Beerwart

Add in info please

[Comment](#)

IN PROGRESS 1

 EQUIPMENT MAINTENANCE EVENT: COOLING SYSTEM OF MX3182

RESOLVE

CUSTOMER: ACME Demo

WORKSITE: Houston , TX

UNIT ID: MX3182

COMP. DESC.: COOLING SYSTEM

ASSIGNED BY : Acme Demo

RECORD MAINTENANCE

[Comment](#)

DONE 1

 EQUIPMENT MAINTENANCE EVENT: TRANS-AUTO OF MX3182

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ALL ACTIONS

+ NUEVA ACCIÓN

Action	Status	Resolution	Assigned To	
 COMPONENT OVERDUE FOR SAMPLING: HYDRAULIC OF 632407	New	Unresolved	Demo Abc	CANCEL REASSIGN ▼
 EQUIPMENT HAS MISSING INFO: ENGINE OF 366215	New	Unresolved	Acme Demo	CANCEL REASSIGN ▼
 EQUIPMENT HAS MISSING INFO: ENGINE OF 632407	New	Unresolved	Acme Demo	CANCEL REASSIGN ▼
 EQUIPMENT MAINTENANCE EVENT: COOLING SYSTEM OF MX3182	In Progress	Unresolved	Acme Demo	REASSIGN ▼
 EQUIPMENT MAINTENANCE EVENT: TRANS- AUTO OF MX3182	Complete	Changed Oil	Acme Demo	REOPEN ▼

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+ Add Filter

ALL ACTIONS

Action	Status	Resolution	Assigned To	
 COMPONENT OVERDUE FOR SAMPLING: HYDRAULIC OF 632407	New	Unresolved	Demo Abc	CANCEL REASSIGN ▼
 EQUIPMENT HAS MISSING INFO: ENGINE OF 366215	New	Unresolved	Acme Demo	CANCEL REASSIGN ▼
 EQUIPMENT HAS MISSING INFO: ENGINE OF 632407	New	Unresolved	Acme Demo	CANCEL REASSIGN ▼
 EQUIPMENT MAINTENANCE EVENT: COOLING SYSTEM OF MX3182	In Progress	Unresolved	Acme Demo	REASSIGN ▼
 EQUIPMENT MAINTENANCE EVENT: TRANS-AUTO OF MX3182	Complete	Changed Oil	Acme Demo	REOPEN ▼

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GESTIONAR ACCIONES

Castrol
LABCHECK
Featuring Labcheck Intelligence

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+ Add Filter

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+ NUEVA ACCIÓN

Action	Status	Resolution	Assigned To	
 EQUIPMENT HAS MISSING INFO: ENGINE OF 366215	New	Unresolved	Demo Abc	CANCEL REASSIGN
 COMPONENT OVERDUE FOR SAMPLING: HYDRAULIC OF 632407	New	Unresolved	Demo Abc	CANCEL REASSIGN
 EQUIPMENT HAS MISSING INFO: ENGINE OF 366215	New	Unresolved	Acme Demo	CANCEL REASSIGN
 EQUIPMENT HAS MISSING INFO: ENGINE OF 632407	New	Unresolved	Acme Demo	CANCEL REASSIGN
 EQUIPMENT MAINTENANCE EVENT: GEARBOX OF MX3182	New	Unresolved	Lori_BPNew Beerwart	CANCEL REASSIGN

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Creado el greater than 12/01/2025 X + Add Filter

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+ NUEVA ACCIÓN

Action	Status	Resolution	Assigned To	
 EQUIPMENT HAS MISSING INFO: ENGINE OF 366215	New	Unresolved	Demo Abc	CANCEL REASSIGN ^
Created By Acme Demo	Customer ACME Demo	Worksite Houston , TX	Unit Id 366215	Comp. Desc. ENGINE
Please add missing information				
Comment				
 COMPONENT OVERDUE FOR SAMPLING: HYDRAULIC OF 632407	New	Unresolved	Demo Abc	CANCEL REASSIGN v
 EQUIPMENT HAS MISSING INFO: ENGINE OF 366215	New	Unresolved	Acme Demo	CANCEL REASSIGN v
 EQUIPMENT HAS MISSING INFO: ENGINE OF 632407	New	Unresolved	Acme Demo	CANCEL REASSIGN v
 EQUIPMENT MAINTENANCE EVENT: GEARBOX OF MX3182	New	Unresolved	Lori_BPNew Beerwart	

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CREATE ACTION(S)

DISMISS ALERT(S)

☐	ALERT TYPE	DETAILS	STATUS	CREATED ON	CUSTOMER	WORKSITE	UNIT ID	COMP. DESC.	
☒		Component Overdue for Sampling: ENGINE of 672547	New	05/02/2023	Acme Demo	Houston, TX	672547	ENGINE	
☐		Component Overdue for Sampling: ENGINE of 631573	New	31/01/2023	Acme Demo	Houston, TX	631573	ENGINE	
☐		Component Overdue for Sampling: ENGINE of 409346	New	31/01/2023	Acme Demo	Houston, TX	409346	ENGINE	

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The tested viscosity is below the range for the oil type specified on the sample information label. Possible causes for viscosity decrease could be the presence of one or more contaminants or mislabeled / improper oil grade. --RECOMMENDATIONS-- Test results indicate one or more tests slightly outside of normal limits. Resample at regular interval.

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Castrol LABCHECK
Featuring Labcheck Intelligence

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RECOMENDACIONES

Mantenimiento: 202509222143
Evaluated by: JHIBBS

The tested viscosity is below the range for the oil type specified on the sample in outside of normal limits. Resample at regular interval.

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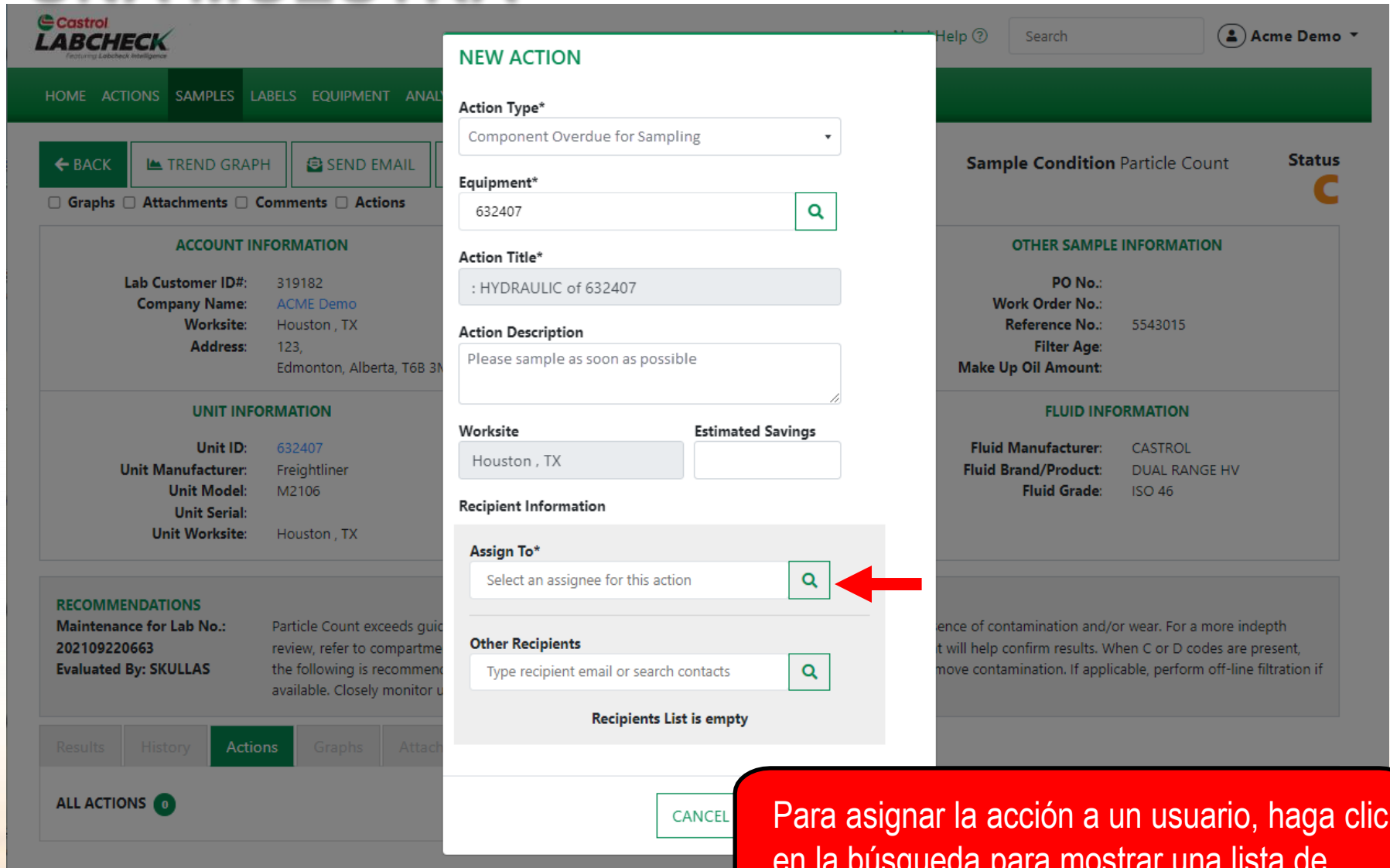
NEW ACTION

Action Type*
Component Overdue for Sampling

Equipment*
632407

Action Title*
: HYDRAULIC of 632407

Action Description
Please sample as soon as possible

Worksite
Houston , TX

Estimated Savings

Recipient Information

Assign To*
Select an assignee for this action

Other Recipients
Type recipient email or search contacts

Recipients List is empty

CANCEL

ACCOUNT INFORMATION
Lab Customer ID#: 319182
Company Name: ACME Demo
Worksite: Houston , TX
Address: 123, Edmonton, Alberta, T6B 3M

UNIT INFORMATION
Unit ID: 632407
Unit Manufacturer: Freightliner
Unit Model: M2106
Unit Serial:
Unit Worksite: Houston , TX

RECOMMENDATIONS
Maintenance for Lab No.: 202109220663
Evaluated By: SKULLAS
Particle Count exceeds guideline. For a more indepth review, refer to compartment the following is recommended. Closely monitor u

OTHER SAMPLE INFORMATION
PO No.:
Work Order No.:
Reference No.: 5543015
Filter Age:
Make Up Oil Amount:

FLUID INFORMATION
Fluid Manufacturer: CASTROL
Fluid Brand/Product: DUAL RANGE HV
Fluid Grade: ISO 46

Sample Condition Particle Count **Status** C

HOME ACTIONS SAMPLES LABELS EQUIPMENT ANALYSIS

← BACK TREND GRAPH SEND EMAIL

Graphs Attachments Comments Actions

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RECOMENDACIONES
Mantenimiento: 202509222143
Evaluado por: JHIBBS
The tested viscosity is below the range for the oil type specified on the sample in outside of normal limits. Resample at regular interval.

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Comentario

Please update the sample as soon as possible

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The tested viscosity is below the range for the oil type specified on the sample information label. Possible causes for viscosity decrease could be the presence of outside of normal limits. Resample at regular interval.

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The tested viscosity is below the range for the oil type specified on the sample information label. Possible causes for viscosity decrease could be the presence of contaminants or oil degradation. Resample at regular interval.

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